

Updating the firmware of a PTC10

December, 2021

To update the firmware in your PTC10 you will first need to email SRS Tech Support (info@thinkSRS.com) with your serial number and a picture of the "About" screen (accessed from the System Setup screen). We will then reply back to you with an email which will has attachments. Copy the attached files to a new USB stick and follow the instructions below.

- 1) Plug the memory stick into the PTC10. After about 15 seconds, a "Firmware update" window will appear on the screen.
- 2) Wait about 10 more minutes for the "Firmware update was successful" message to appear.
- 3) Unplug the memory stick and erase it so that you don't accidentally reprogram the PTC10 again.
- 4) Restart the PTC10.